

Development of a single-tube, 8-color PCR assay for quantifying a *KRAS* mutation panel using true single-molecule counting.

Mike Hoang*, Michael A. Balamotis*, Haeun Lee*, Vladyslava Ratushna*, Andras Saffics*, Mehdi Molaei*, Eleen Y. Shum*, H. Christina Fan*

* Countable Labs, Inc. 1810 Embarcadero Rd. Suite 200 Palo Alto, California 94303

Abstract

KRAS mutations are among the most common drivers of malignant tumors in lung, colorectal, and pancreatic cancers. Identifying specific mutations in solid tumors guides treatment selection, while quantifying mutant-to-wild-type alleles in liquid biopsies helps assess therapeutic response and recurrence risk. Accurate mutation profiling is essential for developing *KRAS*-targeted therapies.

However, standard qPCR assays detect mutations only at ~1% mutant allele frequency (MAF) and require multiple reactions, increasing cost and reducing scalability. Digital PCR (dPCR) improves the sensitivity to ~0.2% MAF but often groups different targets into shared channels, limiting specificity. For low-abundance, limited-input samples such as cell-free DNA (cfDNA), a more sensitive and efficient assay is urgently needed that addresses both sensitivity and specificity challenges.

Here, we present a robust PCR-based method for simultaneous quantification of multiple *KRAS* mutations in a single reaction.

Method, cont.

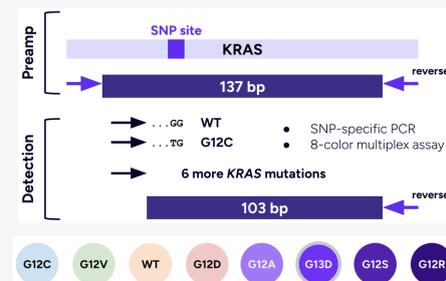


Figure 2. 8-color assay design. Half-nested PCR primer sets were designed to first amplify the *KRAS* locus (preamp) independent of SNP status in a bulk reaction (purple arrows). Once amplified, samples were taken directly into Countable PCR for SNP-specific detection using an 8-color multiplex assay (the black arrows with the purple reverse). The *KRAS* targets are listed by amino acid mutation within a colored circle (optical signature).

A streamlined workflow with the Promega Maxwell® CSC makes cfDNA processing easier that is compatible with Countable PCR.

Total cfDNA was extracted from human blood plasma using a magnetic bead-based automated system (Rapid ccfDNA Kit on the Maxwell® CSC system) to ensure consistent recovery from low-input material.

Single molecule isolation and counting with Countable PCR.

Countable PCR creates ~30 million isolated compartments by centrifugation to enable true single-molecule quantitation. Each target molecule is contained and amplified independently in its own compartment, eliminating amplification bias for multiplex assays, and enabling direct counting.

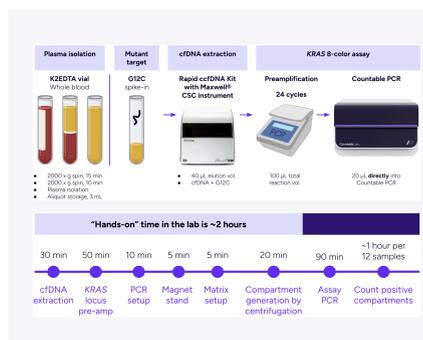


Figure 3. A complete workflow. The workflow begins with plasma isolation from whole blood, synthetic DNA spike-in at a chosen %MAF, cfDNA extraction using the Rapid ccfDNA Kit on the Maxwell® CSC system, *KRAS* target pre-amplification to boost sensitivity, and the *KRAS* 8-plex assay run on Countable PCR.

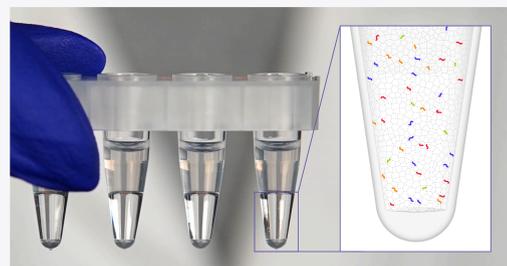


Figure 4. Schematic illustrating the single-molecule isolation of 4 targets into ~30 million picoliter-sized compartments.

Results

Each of the 8-color targets in the assay are detected by a unique optical signature.

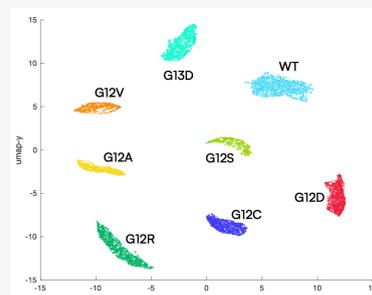


Figure 5. Each target color is distinctly identified. The optical signature in each positive compartment of the Countable matrix is projected onto a 2-dimensional representation using Uniform Manifold Approximation and Projection (the UMAP) analysis. Compartments with similar signatures are placed into the same cluster and are called to a type based on a training set of single-targets that define the map. As shown, each target in the 8-color assay is clearly separated, allowing for determination of counts of each target based on the the number of compartments assigned to a cluster.

Results, cont.

8-color assay achieves the expected performance in counting and specificity against different mutants.

KRAS mutant-specific synthetic DNA templates were tested for on-target specificity as a 1-color or an 8-color assay. To show no counting bias, counts from a 1-color Countable PCR for each mutation should match the counts for that same mutation in the 8-color, and the NTC counts for all targets should be as low as possible relative to real counts.

Table 1. For each mutation tested using a single template in either a 1-color or 8-color PCR reaction, counts are shown per reaction. Along the diagonal (purple boxes), we expect to see counts consistent with the 1-color reaction by *KRAS* mutation type.

1-Plex Counts								
	G12C	G12S	G12V	G12D	G12A	G12R	G13D	WT
G12C	3,097	1,769	1,993	3,363	2,107	3,060	3,446	5,712
8-Plex Counts								
	G12C	G12S	G12V	G12D	G12A	G12R	G13D	WT
G12C	3,149	7	1	0	6	5	0	3
G12S	1	1,792	1	2	0	0	1	19
G12V	0	6	2,059	1	8	1	0	4
G12D	0	2	5	3,701	13	0	0	18
G12A	0	1	4	1	2,115	1	1	1
G12R	1	1	0	0	0	3,922	1	5
G13D	0	1	0	2	0	1	4,016	22
WT	4	4	1	4	0	2	105	5,325
NTC Counts								
	G12C	G12S	G12V	G12D	G12A	G12R	G13D	WT
	12	3	3	3	21	5	7	10

Determining sensitivity of *KRAS* detection in extracted samples.

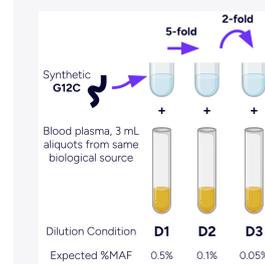


Figure 6. Dilution scheme. Total cfDNA was extracted from 3 mL of human plasma using the Rapid ccfDNA Kit on the Promega Maxwell® CSC system. Aliquots were spiked-in with dilutions from a synthetic *KRAS* G12C template to create an estimated %MAF over three points of diminishing abundance, starting from an expected MAF of 0.5% in sample D1.

Countable PCR confidently detects 0.08% MAF for G12C in cfDNA.

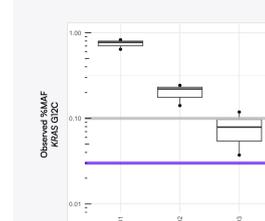


Figure 7. Pre-amplification of cfDNA boosted the number of mutant targets above background. Support of a broad counting range allowed the Countable PCR *KRAS* assay to detect rare targets better than existing methods (purple vs. gray LoD line).

The broad counting range of Countable PCR allowed the detection of >140,000 WT *KRAS* molecules (Figure 7 and Table 2) against ~113 G12C counts in a single reaction, an average MAF of 0.08% and above the experimentally determined limit of detection (purple line, an average of 44 counts, N=8), in contrast to dPCR assays (gray line) that cannot reliably detect below 0.1% MAF.

Table 2. Molecule counts per reaction for each dilution, shown as average value of technical replicates. Average counts are for wild-type and G12C mutant after pre-amplification from 25 ng of total cfDNA. The expected copy input of G12C was ~5 copies per reaction into pre-amplification.

Sample	Dilution	WT		G12C		%CV	Replicates
		Ave. counts	Ave. counts	Ave. %MAF	%CV		
D1	1-fold	143,030	1,072	0.75%	12.8%	N=3	
D2	5-fold	109,247	216	0.20%	26.5%	N=3	
D3	10-fold	147,206	113	0.08%	51.9%	N=3	
Blank	-	-	18	-	-	N=8	

Conclusion

The Countable PCR single tube *KRAS* 8-color assay can detect 0.08% MAF for the G12C mutation in human cfDNA samples.

We unambiguously identified seven *KRAS* mutants in a single reaction. When coupled with a pre-amplification step, the assay showed detection sensitivity to 0.08% for the G12C mutation, approximately one order of magnitude more sensitive than published assays. Results were reproducible across technical replicates and were achieved with minimal optimization.

We developed an 8-color assay for *KRAS* mutations and wild-type, all in one reaction

Using pre-amplification to boost target counts, we measured over 140,000 targets in the same reaction, demonstrating how a broad dynamic counting range supports sensitive target detection

Countable PCR is compatible with the Rapid ccfDNA Kit on the Maxwell® CSC system for cfDNA extraction

Using *KRAS* cancer mutations as an example, we showed that Countable PCR enables profiling of multiple targets in a single experiment with high sensitivity and reduced time for optimization. With straightforward and sensitive 8-color multiplexing, this approach opens new avenues for research and diagnostic panels targeting other cancers.

Background

99% of known *KRAS* mutations are covered by 7 exclusive single base pair mutations

Some *KRAS* mutations can be drugged (G12C, G12V, G12D), while other mutations have yet to be treatable

Knowing a specific *KRAS* mutation helps guide treatment options

Countable PCR overcomes the multiplexing limitations of qPCR and dPCR, enabling precise quantification of gene panels in a single reaction, providing a practical solution for biomarker validation post NGS studies.

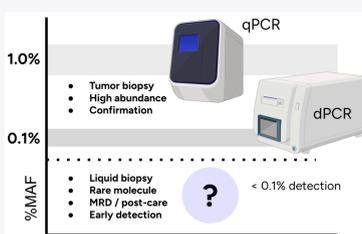


Figure 1. Rare target discovery is limited by detection sensitivity. *KRAS* assays that rely on qPCR and dPCR have limited sensitivity, reducing their effectiveness in liquid biopsy (cfDNA) screening, post-treatment minimal residual disease (MRD) screening, and early detection testing.

Method

Countable PCR 8-color *KRAS* assay uses a pre-amplification step to improve the sensitivity of detection.

We utilized Countable PCR, which enables direct counting of single molecules across a broad dynamic range, to develop an 8-color assay that differentiates and quantifies both common and rare *KRAS* mutations in a single reaction, with higher precision and sensitivity than other platforms. The inclusion of a SNP-agnostic pre-amplification step accurately quantifies abundant wild-type and low-frequency mutant allele without loss of accuracy. Each *KRAS* mutation (G12C, G12V, G13D, G12D, G12A, G12S, G12R), along with the wild-type allele, is represented by one of eight unique optical signatures.

